LED-Lampenhaus
LED Lamp House

MLH2-Series

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Vielen Dank

General Information

This high quality product has been designed and developed to require a minimum of maintenance. Please follow the guidelines below and read this manual in its entirety to ensure safe operation of the unit.

Symbols

The following symbols are used in this manual.

CAUTION! This symbol indicates a potential risk and alerts you to proceed with caution.

CAUTION! This symbol indicates a potential danger of optical radiation.

Warning

- Please always disconnect mains plug before cleaning the device.
- For indoor use only.

Safety Information

- To avoid eye damage, don't look directly into the light source when turned on.
- Versions with UV light release intensive ultraviolet radiation during operation.

A major part of the light is invisible. The beams can damage your health. Protect your eyes and skin during operation.

Important: The evaluation of eye safety occurs according to the standard CIE S009/E:2002 („photo biological safety of lamps and lamp systems“). Within the risk grouping system of this CIE standard, the LEDs specified in this data sheet fall into the „low risk“ group (relating to devices in the visible spectrum with an exposure time of 100s). Under real circumstances (for exposure time, eye pupils, observation distance), it can be assumed that no endangerment to the eye exists from these devices. As a matter of principle, however, it should be mentioned that intense light sources have a high secondary exposure potential due to their blinding effect. As is also true when viewing other bright light sources (e.g. headlights), temporary reduction in visual acuity and afterimages can occur, leading to irritation, annoyance, visual impairment, and even accidents, depending on the situation.

Function and Operation

Package Contents

LED Lamp House MLH2	1
LED Brightness Control Unit (TC)	1
Multi Adapter Power Supply	1
User Manual	1

Technical Data

Luminous flux	max. 1.250 lm (PW)
Illuminant	30 W Power-LED
LED life span	50.000 h (according to manufacturer)
Power consumption	approx. 36 W
Input power supply	100 - 240 V AC, 50 - 60 Hz
Output power supply	24 V DC, 1 A
Dimensions (L×W×H)	100 mm × 84 mm × 100 mm
Material	anodized aluminium
Color	titanium and black
Weight	approx. 1,5 kg
Ambient temperature	10°C to 35°C
Cooling system	passive cooling without fan

Versions

MLH2-PW pure white (5.700 K), CRI 70	100-011163
MLH2-NW natural white (4.000 K), CRI 80	100-010945
MLH2-WW warm white (3.000 K), CRI 80	100-011176
MLH2-PC pure white (5.700 K), CRI 90	100-011174

Accessories

Adapter for Zeiss Microscopes	100-011008
Adapter for Leica Microscopes	100-011030
Adapter for Olympus Microscopes	100-011183
Adapter for Motic Microscopes	100-011010
Adapter for Nikon Microscopes	100-011012
Adapter for Nikon Microscopes (bayonet)	100-011180

Getting Started

The LED lamp is a plug & play device very easy to use.


Please check before use whether...

...the mains plug is connected.

...the device is switched on (the blue LED light glows).

...the brightness is set to the desired value.

Description of Buttons and Switches

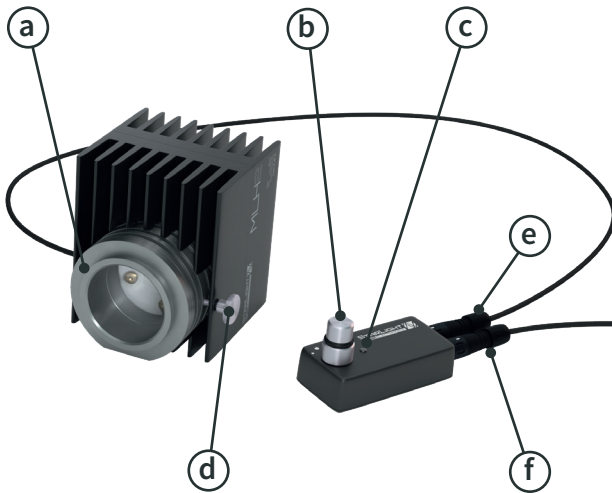- a) **Adaptor**
To mount on microscope
- b) **On/Off Switch and Brightness Control**
Activation of the memory function (see chapter below)
- c) **Operation Control**
The blue LED glows as soon as the ring light has been switched on
- d) **Knurled Screw**
To adjust the focus distance (0 - 12 mm)
- e) **Connection from the LED Lamp House**
Please connect correctly before use
- f) **Connection from the power supply**
Please connect correctly before use

Brightness Control with Memory Function

The brightness can be processor controlled adjusted. The highest level is indicated by a double blinking of the LED.

Memory function:

1. When switching off the device by pushing the button the last setting is automatically saved for the next use.
2. Before an interruption of the power supply (e. g. withdrawing of the power adaptor or switch off of the main switch) the last setting can be saved individually by the user:
 - At the preferred level please push the button for 3 seconds.
 - A double blinking of the LED signals a successful storing.
 - Now, the power supply can be interrupted.

Technical Assistance

Troubleshooting

The information in the table below can help to resolve the possible operating problems quickly.

Problem	Causes	Solution
No function at the device	no network connection	Please check whether the power adaptor is connected.
	electronic defective	Please contact your distributor.
	LED defective	Please contact your distributor.
LED gets dark	thermal overload of the system	Please ensure an ambient temperature from 10° to 35°C.
	LED life span exceeded	Please contact your distributor.

Storage, Maintenance, Repair

The device is maintenance-free. Some points, however, you should take into consideration.

1. Please unplug the device before cleaning.
2. For the cleaning of the housing please use a smooth, semi-wet cloth, if necessary with a smooth household cleaning fluid.
3. No moisture should entering into the housing interior.
4. To avoid damages during storage or transport please retain the packaging.

Repairs are only allowed by trained and experienced personnel. In case of improper handling the warranty expires.

Technical Service

Web Resources

Visit our StarLight website at starlight-sl.com for:

- Contact details of technical customer service.
- StarLight product overview and price list as well as infos about accessories and spare parts.
- Additional product information and special offers.

Contact

For information or technical assistance contact your local distributor or directly StarLight Opto-Electronics GmbH & Co. KG.

Warranty

StarLight Opto-Electronics GmbH & Co. KG warrants that this product will be free from defects in material and workmanship for a period of one (1) year from date of purchase. If a defect is present, StarLight will, at its option, repair, replace, or refund the purchase price of this product at no charge to you. Precondition is the device is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products.

It is expressly agreed that this warranty will be in lieu of all warranties of fitness and in lieu of the warranty of merchantability.

Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead, it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you