


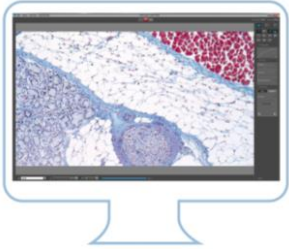

Stability –
giving you a reliable research tool
you can count on.



UPDATE INFORMATION - PROGRES GRYPHAX® microscope camera software

We are glad to provide you with our new PROGRES GRYPHAX® software version 1.1.8. for MAC OS. It is free of charge. You can download the latest software version from download page below after registration.

JENOPTIK PROGRES GRYPHAX® microscope camera software

[DOWNLOAD UPDATE](#)

SHORT LIST OF NEW FEATURES

- New feature "Tree View" to save, organize and handle media destination folder structure
- Z-Stacking - New option to change combining mode (fast or parallax correction for stereo microscopes)
- New option for fast image record, to shorten image saving time
- New Video storage format "mp4" and "webm" for Mac OS
- Multi-fluorescence tool available during GUI scaling
- Added option to change zoom factor for magnifier tool
- Added option to change size of Gallery thumbnails
- New storage option "PNG 16 bit"
- Gallery and "Side by Side" view revised and updated
- Added software "Service Mode" for trouble shooting
- Added new [USB 3.0 Camera Service Tool](#) for GRYPHAX® cameras to update firmware and add hot pixel correction

SHORT LIST OF ADDED FEATURES from previous version

- Added camera support for PROKYON, KAPELLA and RIGEL
- New optimized work flow for Multi-Fluorescence tool
- New option to password protect "User profiles"
- New stereo microscope shift compensation for Z-Stacking tool
- Enhanced performance for Panorama tool
- Added higher software scaling factors for high resolution displays (e.g. 4k/ 5k resolution)
- Added new image resolution [3.7 Mpixel](#) (2560 x 1440 px) for ARKTUR cameras
- Added measurement for panorama, z-stacking and fluorescence images from gallery
- Added support for export of image Meta-data and software settings
- New feature "[Ocular view](#)" to adapt the color impression from ocular to images on screen

BUG FIXES & IMPROVEMENTS

- Fixed an issue for Panorama image stitching
- Fixed an issue for scale bar lengths calculation
- Measurements description lengths reduced
- Fixed an issue that Print was not working on certain computer
- Fixed an issue for colour reproduction of NAOS camera (Firmware update by service tool necessary)
- Fixed an issue that caused warning after software download from internet
- Fixed an issue that recorded image where shown rotated on certain operating systems
- Fixed an issue for image analyze & send file option

PREVIEW for upcoming versions

- Update of Measurements

PROGRES GRYPHAX® software version 1.1.8 is available for the following platforms and operating systems

- macOS Sierra (10.12) - 64 bit
- macOS El Capitan (10.11) - 64 bit

Important note:

- GRYPHAX Mac version 1.1.8 is the last software version with ProgRes® Firewire camera support.

It is supporting the following ProgRes® and PROGRES GRYPHAX® cameras

- PROGRES GRYPHAX® PROKYON
- PROGRES GRYPHAX® KAPELLA
- PROGRES GRYPHAX® RIGEL
- PROGRES GRYPHAX® NAOS
- PROGRES GRYPHAX® ARKTUR
- PROGRES GRYPHAX® SUBRA
- ProgRes® C14^{Plus}
- ProgRes® CF^{cool}
- ProgRes® MF^{cool}

[LEARN MORE](#)

[VIDEO TUTORIALS](#)

[HELPFUL? WE APPRECIATE YOUR FEEDBACK](#)

[FEEDBACK](#)

We looking forward to hearing from you and remain sincerely,

Your PROGRES GRYPHAX® Team