




User Guide for PROGRES GRYPHAX® MetaMorph driver

The “MetaMorph driver” for PROGRES GRYPHAX® USB 3.0 cameras enables user to operate with MetaMorph microscopy automation & image analysis software solutions from Molecular Devices, LLC. The driver delivers essential functionalities of PROGRES GRYPHAX® cameras. Driver is supported under Windows operating systems.

The GRYPHAX MetaMorph driver is part of PROGRES GRYPHAX® software installation (as of version 2.0.0 or newer).

General Preparation:

The MetaMorph software version **7.8.0.0 – 32 bit** has to be installed before installation of GRYPHAX MetaMorph driver. Recommended USB 3.0 interface card (with Renesas chipset) has to be installed in advance!

Overview:

GRYPHAX camera operating under MetaMorph driver:

