




**Stability –**  
giving you a reliable research tool  
you can count on.



## User Guide for Report tool of PROGRES GRYPHAX® software

The “[Report tool](#)” of PROGRES GRYPHAX® software enables user to review and export measurement data, software settings and image meta-data directly at GRYPHAX software.

### General description:

The “[Report tool](#)” is part of the GRYPHAX software. It enables user to review and export measurement data from live and recorded images, current software settings and media files meta-data directly at GRYPHAX software.

### Preparation:

To use all feature of *Report tool*/you have to calibrate your microscope objectives in advance! Otherwise the *Report tool* cannot be display measurement data information. To create measurement calibration you have to open GRYPHAX Preferences and navigate to option “Device Configuration”.

To calibrate the microscope you can use the PROGRES GRYPHAX® [calibration slide](#) (order number: 563683).



### Overview:

Report tool contains the data tables and views for:

- Measurement data & statistics
- Image meta-data view
- Software settings view

Show image information from selected media file

